FFC-100 Fiber Frequency Comb Quick Start Guide

Model No. FFC-100-100 and FFC-100-200 Document Last Updated on 2023/06/08 22:32

Please read Limited Warranty and General Warnings and Cautions prior to operating the FFC-100.





FFC-100 Fiber Frequency Comb

Useful Links

Click here for the Vescent manuals page. Click here for the FFC-100 API. Click here for the SLICE-FPGA-II Locking Manual Click here for the FFC-100 web page. Click here for the Github page for FFC-100 firmware revisions Click here for the FFC-100 Manual Please check back for added functionality. Contact sales [at] vescent [dot] com for questions and corrections, or to request added functionality.

Notice



Do not block the airflow vents on the side of the chassis or the fan inputs & outputs on either the FFC-100 or the SLICE-FPGA.

The mode-lock indicator on the GUI front panel cannot detect CW breakthrough. Keep the oscillator current within the range specified in your product's final test documentation or CoC.

Operating the FFC

This document provides instructions on how to operate the Vescent Photonics FFC-100 when controlled by the Vescent SLICE-FPGA or SLICE-FPGA-II.

Purchase Includes

- FFC-100 rack-mountable Fiber Frequency Comb
- Power cord for your country (if known)
- Certificate of Conformance

Absolute Maximum Ratings and Power Input

Note: Reported specs are under laboratory conditions. Performance degradation can occur in deployed operation depending on the environment

Parameter	Rating
Environmental Temperature	15°C <t<30°c< td=""></t<30°c<>
Environmental Humidity	<60%
Environmental Dew Points	<15°C
Maximum AC Line Input Current	2 A
Tab. 1: Absolute Maximum Ratings	· · · · ·

The FFC-100 employs a proprietary design hybrid power supply that is both low noise and capable of accepting a range of AC input line voltages. It will accept input line voltages within the ranges shown in table 2.

Parameter	Value	Units
Input Line Voltage	100-240	VAC
Frequency	50-60	Hz
Phase	1 phase	-
User-serviceable fuse ¹⁾	T 2.0 A L 25	50V
Tab. 2: Input Voltage Specific	ations	

Proper Usage



If this instrument is used in a manner not specified by the manufacturer in this manual or other relevant literature, protection provided by the instrument may be impaired.

Successful implementation of the FFC-100 depends critically on the design of the whole system: FFC-100, phase locking electronics, and any references to which the FFC is locked or *vice versa*.

Initial Set-up

- 1. Ensure all cables and fibers are connected according to the connection diagram figure 2.²⁾
- 2. Ensure the RF input amplitudes to the FPGA are between -30 dBm and -15 dBm. If power levels are higher than -15 dBm, utilize an appropriate RF attenuator.

- 3. Turn on the power to the Fiber Frequency Comb (FFC) and the FPGA, as well as the reference laser.
- 4. Ensure the optical input power to the fOPT input port is < 1.5 mW.





Parameter	Rating
f _{opt} Power Input	<1.5 mW max
Tab. 3: Operating Parameters	

FFC-100 Operation

FFC Start Up GUI

1. Touch the "System Off" button at the top of the front panel screen (figure 3) to bring up a dropdown menu. Select "Standby" (figure 4) to turn on temperature control of the pump diodes and oscillator cavity.

<u>*</u>	SYSTEM OFF	Mode-locked		
	OSC CAV	ITY		
	PZT Voltage [V]: 0.00 0 DC Bias IVI	100 Setpoint I°C1 Actual I°C1		
	PZT 0.00 Temperate	ure 0.000 0.000		
	PUMP LAS	SERS		
	OSC Mod. [mA]: 0 -100	100	1	
	Current [mA] Temp ["C]	Current [mA] Temp [°C]		
0	OSC 0 0.000 A	AMP 0 0.000	Fig. 3: The FFC-10	0 GUI in off state
	🗶 🗖 sys			
	STA			
			100	
		ER ON	Setpoint ["C] Actual ["C]	
		0.00 Temperature	0.000 0.000	
	Ů	PUMP LASERS		
	OSC Mod. [m/	A]: 0 -100	100	
		rent [mA] Temp [*C]	Current [mA] Temp [*C]	
		0 0.000 AMP	0 0.000	

Fig. 4: Select Standby laser state

 Once the temperature servos have stabilized and "Standby" has stopped blinking, touch the "Standby" drop-down menu (figure 5) and select "Laser On" (figure 6). The FFC should commence lasing and the Mode-locked indicator should turn green (figure 7).



Fig. 5: The FFC-100 GUI: in Standby

mode

2	SYSTEM OFF	Mode-locked
	STANDBY	CAVITY
9	PZT	100
5	LASER ON	Setpoint ["C] Actual ["C]
	PZT 21.00 Temp	erature 25.000 0.000
	PUM	P LASERS
	OSC Mod. [mA]: 0 -100	100
	Current [mA] Temp ["C]	Current [mA] Temp [°C]
	OSC 50 0.000	AMP 70 0.000
Fig. 6: The FFC-100 GUI: S	Standby to System On	
2	LASER ON	Mode-locked
	osc	CAVITY
	PZT Voltage [V]: 21.00 0	100
	DC Bias [V]	Setpoint [°C] Actual [°C]
	P21 21.00 Temp	erature 25.000 0.000
	PUM	P LASERS
	OSC Mod. [mA]: 0 -100	100
	Current [mA] Temp ["C]	Current [mA] Temp [°C]

Fig. 7: System On

 If necessary, adjust oscillator current by either holding down the "Osc" touch screen button and typing in a desired set-point (figure ##), or briefly pressing the button and using the two control knobs (left changes digits, right adjusts values) to set a desired value.



Fig. ##: Adjust oscillator current via keypad

- 4. Cavity temperature can be adjusted in the same way.
- 5. Touch the "PZT" touch screen button (figure 8) to bring up a drop-down menu and select "Full Range x20" to enable the PZT servo.





FPGA Control

For locking the FFC-100 with the Vescent SLICE-FPGA-II Module, click here. Otherwise, if using the SLICE-FPGA, follow the instructions below.

The SLICE-FPGA dual-channel Offset Phase Lock Servo can be used to phase lock f_{CEO} to a reference (TBD) and f_{opt} to a reference laser such as the Rio Planex.

If you have not already done so, install Python and the FPGA software for controlling the SLICE-FPGA dual Offset Phase Lock Servo.

Software Startup

- 1. Open a WinPython command window and navigate to the folder "GUI and Firmware".
- 2. Start the SLICE-FPGA control GUI by typing "python XEM_GUI3_VPv4.py", figure ##).



Fig. ##: Start FPGA GUI from Terminal

3. A start-up menu should appear (figure ##). Make sure "superlaserland_v12.bit" is selected and

select an appropriate clock option. Press OK.

C:\WINDOWS\system32\cmd.exe - python XEM_GUI3_VPv4.py						
C:\Users\Bennett\Documents\NIST FPGA Software\Wi \NIST FPGA Software\GUI and Firmware	inPython-64bit-2.	7.10.3\python-2.7.10.amd64>cd C:\	Users\	Bennet	t\Documer	nts ^
C:\Users\Bennett\Documents\NIST FPGA Software\GU XEM_GUI3 - INFO - STAR XEM_GUI3.SLL_JD2 - INFO - Inst Http server launched Starting httpd server Ø {0: '1624000EUI'} 1624000EUI	UI and Firmware≻p RTING MAIN FUNCTI tantiating Superl	oython XEM_GUI3_VPv4.py CON aserLand class				
10	Initial configuration		?	×		
Con Firm ()	nnected FPGAs mware file Send firmware Internal clock (100 MHz)	Name = , Serial # = 1624000EUI, Color = superlaserland_v12.bit O Connect to an already running box (NOT REALLY W External clock (200 MHz, divided by 2 internally)	ORKENG Y	• • ET)		
		OK	Cancel			
						Ŷ

Fig. ##: FPGA GUI Start Screen

4. Navigate to the "Filter Settings" tab and select "Narrowband (6MHz)" for both DDC0 and DDC1 (figure ##).

۰.				
-	CED Lock OpricaliLock Counters Differ Se DIC settings	the settings	Filter Settings	
		Narr	owband (6 MHz)	
		\sim		
	COC 9 New BW.		Iserconbend (5 Mez)	O Webard PRI (S0 MH3
	COC 1 film BH	 Wideband (25 M-Ir) 	Harronbard (6 Nict)	O Welcard Fill (SD HH)

Fig. ##: Set Loop Filter Bandwidth

Locking f(CEO)

 Navigate to the "CEO Lock" tab and adjust the "Offset DAC 0" slider near the top left by clicking and dragging the slider until the the data in the Baseband IQ plot is circular (figure ##). This centers the beat note near the reference frequency. It is also possible to adjust the FFC-100 oscillator current on its front panel to make this adjustment.



Fig. ##: Centering f(CEO) on reference frequency (Baseband IQ optimization)

2. Press the "Lock" button (figure ##, top middle). If the system doesn't lock, change the VCO sign to the opposite polarity (top right) and try again. If the system still won't lock, try lowering the K_p value (bottom left).



Fig. ##: Locking f(CEO)

Locking f(opt)

 Navigate to the "Optical Lock" window. Center the beat note near the reference frequency: adjust the "Offset DAC 1" slider (or your reference laser frequency) until you see a circular Baseband IQ diagram (figure ##).



Fig. ##: Centerig f(opt) on reference frequency (Baseband IQ optimization)

2. Press the "Lock" button (figure ##). If f_{opt} doesn't lock, change the VCO sign to the opposite polarity and try again. If the system continues to not lock, lower the K_p value.\\



Fig. ##: Locking f(opt)

3. Adjust PID settings (bottom middle of figure ##) accordingly to lower the integrated phase noise of each parameter (f_{CEO} and f_{opt}). The default settings provided in the software are a good place to start but tweaking the values can often improve performance.

Slow Loop Feedback

Firmware Version FL-V1.2

Slow loop functionality will be provided in a near-term software update.

- 1. Open your computer's device manager from the control panel and find the COM port number for "STMicroelectronics". Never assume that this number is the same as previous times. (If it doesn't appear, make sure your computer is connected to the FFC.)
- 2. Navigate to the "Slow Loop" tab in the GUI and enter the COM port number into "FFC COM Port".
- 3. The PZT setpoint is set to an intermediate value (50V) to keep the servo from railing. This can be adjusted if desired.
- 4. Click "Activate Temperature Slow Loop".

Firmware Version G2-V1.2

Slow servo functionality is designed to allow indefinite F_{OPT} locking times despite long-term drifts in the reference laser's frequency outside the dynamic range of the FFC's PZT feedback. This is achieved by slow temperature feedback to the oscillator of the FFC, which locks the PZT output voltage to a setpoint voltage near the middle of its dynamic range. To enable the slow servo, first lock FOPT normally, but set the DC offset as close to 3V as possible (PZT output range spans 0-60 V with 10V/V transfer function). Next, navigate to the home screen of the FFC-100 and press the "Slow Servo" button at the top of the screen

	SYSTEM OFF	Slow Servo	Mode-loc	ked 🔵
	PZT Voltage [V]: -1.0	0.0		100.0
5	DC Bias [V]		Setpoint [°C]	Actual [°C]
	PZT 0.00	Temperature	25.000	24.998
Ľ		PUMP LASERS		
	OSC Mod. [mA]: 0.0	-100.0		100.0
	Current [mA]	Temp [°C]	Current [mA]	Temp [°C]
O	OSC 0	0.000 AMP	0	0.000

Fig. 9: The slow servo button is located on the home screen to the right of the System Off/Laser On button.

This will activate the slow servo but does not turn on the feedback. You can tell that the slow servo has been activated by the blue background on the "Setpoint [°C]" button.

S	SYSTEM OFF	Slow Servo	Mode-loc	ked 🔿
Ĥ		OSC CAVITY		
	PZT Voltage [V]: 1.0 0.0			100.0
5	DC Bias [V]		Setpoint [°C]	Actual [°C]
Ā	PZT 0.00 T	emperature	25.000	25.003
Ľ		PUMP LASERS		
	OSC Mod. [mA]: 0.0 -100	.0		100.0
	Current [mA] Tem;	p [°C]	Current [mA]	Temp [°C]
\bigcirc	OSC 0 0.0	DOO AMP	0	0.000

Fig. 10: When the slow servo is activated, the background of "Setpoint [°C]" turns blue

Next, press and hold the "Setpoint [°C]" button until the slow servo menu comes up. The menu allows a user to set the temperature setpoint as normal, and to set the voltage setpoint that the slow servo will try to keep constant by changing the temperature of the oscillator. We recommend setting this to 30V, which is right in the middle of the FFC-100's range. The gain comes preconfigured for each unit, but can also be changed through this menu (WE DO NOT ADVISE changing the gain without input from the Vescent technical staff). The polarity button will toggle the direction that the slow servo adjusts the temperature to keep the setpoint voltage constant (this will be preset by Vescent staff and should not be changed). Setting this to the incorrect value can result in temperature runaway. To change a value, simply touch the field to edit with the rotary knobs, or press and hold to bring up the keypad. To turn the slow servo on, press the "OFF" button. It should change to a white background which says "ON". To turn it back off, simply press the button again.



Fig. 11: The slow servo menu allows the user to change the setpoint, gain, and polarity of the servo, and also to turn it on or off.

While the slow servo is on, the color of the "Setpoint [°C]" button will indicate the difference between the setpoint voltage, and the actual voltage. If the difference is $\leq \pm 6V$, the background will turn green. If it is $\geq \pm 6V$, it will turn yellow, and if it is $\geq \pm 54V$ it will turn red. If the background is blue, this means the slow servo has turned off.

	STANDBY	Slow Servo	Mode-loc	ked 🔵
٦		OSC CAVITY		
	PZT Voltage [V]: -1.0	0.0		100.0
5	DC Bias [V]		Setpoint [°C]	Actual [°C]
Ā	PZT 0.00	Temperature	25.000	24.983
		PUMP LASERS		
	OSC Mod. [mA]: 0.0	100.0		100.0
	Current [mA]	Temp [°C]	Current [mA]	Temp [°C]
0	OSC 0	0.000 AMP	0	0.000

Fig. ##: The slow servo indicating the Voltage is within $\pm 6V$ of the setpoint.

	STANDBY	Slow Servo	Mode-loc	ked 🔵
				100.0
5	DC Bias [V]	.0	Setpoint [°C]	Actual [°C]
$\mathbf{\tilde{h}}$	PZT 0.00		25.000	24.992
	OSC Mod. [mA]: 0.0 -1	PUMP LASERS	1	100.0
	Current [mA] T	emp [°C]	Current [mA]	Temp ["C]
O	OSC 0	0.000 AMP	0	0.000

Fig. ##: The slow servo indicating the Voltage is more than $\pm 6V$ away from the setpoint, but less than $\pm 54V$.

	STANDBY	Slow Servo	Mode-loc	ked 🔿
		OSC CAVITY		
\leq	PZT Voltage [V]: -1.0	0.0		100.0
5	DC Bias [V]		Setpoint [°C]	Actual [°C]
Ā	PZT 0.00	Temperature	25.000	24.982
Ľ		PUMP LASERS		
	OSC Mod. [mA]: 0.0	-100.0		100.0
	Current [mA]	Temp [°C]	Current [mA]	Temp [°C]
0	OSC 0	0.000 AMP	0	0.000

Fig. ##: The slow servo indicating the Voltage is more than ±54V away from the setpoint. This scenario is very unlikely.

Remember to shut off the slow servo, either by disabling it from the home menu or by pressing the On/Off button in the slow servo menu, when unlocking the FFC-100.

1)

Located in power receptacle on rear panel

2)

Single frequency reference laser: for instance a RIO Planex[™] 1550 nm laser.

